

technoclone



CEVERON[®] MFU 500

For Research Use Only


Microparticle filtration unit






diapharma[®]
From Hemostasis to Cell Death

A clear view into the future!

Ceveron[®] MFU 500
enables easy, quick and standardised
removal of *MICROPARTICLES* from
platelet poor plasma.

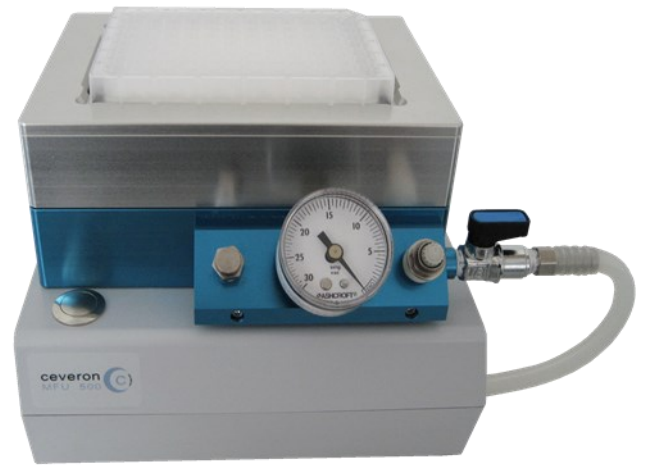

EASY & QUICK

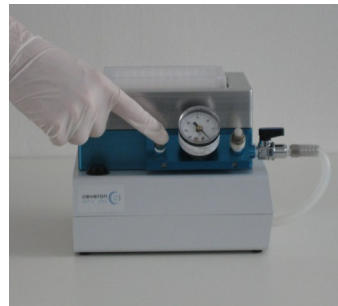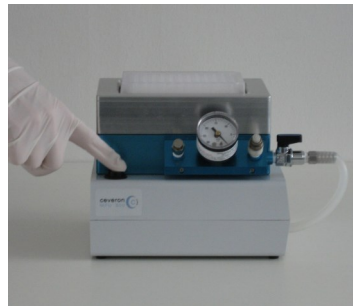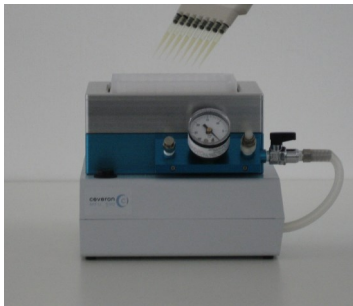

Pipette sample into
the filter plate

Push the "START"
button and filtration
starts automatically

After filtration is
terminated push the
"VENTILATION" button

Filtered sample can be
collected from the
receiver plate


SMALL & COMPACT



Filtration is as effective as
ultra-centrifugation to re-
move thrombogenic MP's
from PPP.

Dimensions MFU 500:
Length: 26.00 cm
Width: 20.00 cm
Height: 16.05 cm
Weight: 3.58 kg

STANDARDIZED

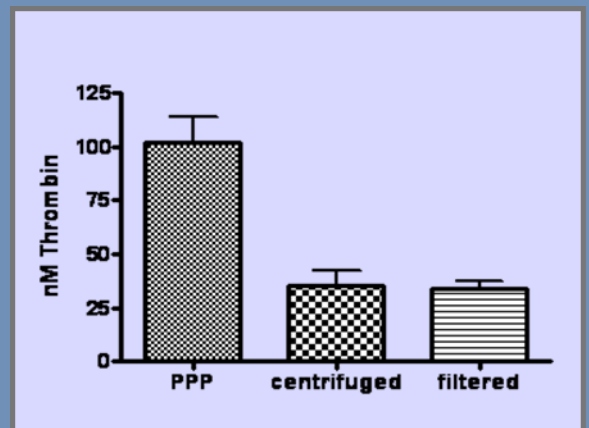

REF 9821010 Ceveron[®] MFU 500 **For Research Use Only**

CEV002E.002 06/2013

DiaPharma Group, Inc.
8948 Beckett Road, West Chester, OH 45069 | USA
Tel: 1-800-526-5224 | Fax: 513-860-9635
info@diapharma.com | www.diapharma.com
Technoclone GmbH
Brunner Str. 67 | 1230 Vienna | Austria
Tel: +43 1 86 373-0 | Fax: +43 1 86373-44
sales@technoclone.com | www.technoclone.com

