

Ecarin Diluent - 20 mL

**ED0413-20****2-8 °C, Do not freeze****Ready for Use**

For Research Use Only. Not for use in diagnostic procedures.

Properties

A liquid stock solution of Tris buffer containing bulking agents for stabilization of Ecarin.
Ready for use.

Composition

0.05 M Tris pH 7.6 (at 25°C), with NaCl, Bovine Serum Albumin (BSA) and polyethylene glycol.

Stability

A sealed vial is stable at 2-8°C until the expiry date printed on the label.

An opened vial is stable for one month at 2-8°C, provided that any microbial contamination is avoided during use.

Application

Buffer for reconstitution and dilution of the purified snake venom prothrombin activation enzyme Ecarin and of the crude venom of Echis Carinatus.

Reconstitute your Ecarin in ED0413-20 Ecarin Diluent according to the package insert for the Ecarin to be used. Use ED0413-20 Ecarin Diluent to further dilute the reconstituted Ecarin to obtain your desired concentration.

Package

Working Solution, 6 x 20 mL



Rossix AB
Taljegårdsgatan 3B
S-431 53 Mölndal, Sweden

www.rossix.com
info@rossix.com

Revision: September 2021