**REF TC41067 PAI-1 Protein ~500 IU****REF TC41069 PAI-1 Protein ~250 IU**

3050100 Rev. 006 [03] 10/2019

**RUO****PAI-1 Protein - English****SUMMARY**

Activity content is determined by PAI-1 Actibind® ELISA Kit, which is standardized against the international PAI-1 standard. Briefly the Actibind assay measures the complex formation between the immobilized t-PA and active PAI-1, detecting complexed PAI-1 with a specific peroxidase labelled antibody.

**REAGENTS**

Plasminogen activator inhibitor type-1 (PAI-1), lyophilized is produced as recombinant protein in *E.coli* (1). The product is lyophilized from 50 mM Sodium acetate, 100 mM sodium chloride, 60 mM L-Arginine-monohydrochloride and 0.01 % Tween 80.

The exact lot specific PAI-1 activity is provided on the vial label.

**Materials required** (not supplied with the product)

- Distilled water
- Test Kit
- **REF** TC16075 PAI-1 Actibind® ELISA Kit

**Warning and precautions**

- For research use only
- Not for drug use
- A Material Safety Data Sheet for this product is available from [www.technoclone.com](http://www.technoclone.com)

**Stability and storage**

The expiry date printed on the labels applies to storage of the unopened bottles at +2...8 °C.

Stability after reconstitution:

≤ -70 °C
2 years or until
expiry date

The reconstituted product should be stored in aliquots.

Repeated freeze thawing is **NOT** recommended.

**TEST PROCEDURE****Preparation of reagent**

- The distilled water and reagent have to reach room temperature (18...25 °C) before mixing.
- The lyophilized product is reconstituted by addition of 1.0 mL distilled water and gentle agitation for 5 minutes at room temperature
- Centrifuge 3 minutes at 1000 g
- Whenever full activity is required, a measure to stabilise its activity has to be taken e.g. addition of plasma, vitronectin, low pH or dibasic aminoacids.

**LIMITATION OF THE TEST**

- Use only clean pipettes.
- Use reconstitution liquid (distilled water) at room temperature. Low temperature might cause flocculation.

**STANDARDISATION**

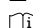
Each lot is traceable to the International Standard for Plasminogen Activator Inhibitor-1 (PAI-1) Plasma.

**LITERATURE**

Please contact Technoclone or your local distributor.

- (1) Ehrlich H.J. et al: Alteration of serpin speciality by a protein cofactor. J. Biol. Chem. 265, 13029-13035, 1990.

**SYMBOL KEYS**

	Manufacturer	<b>RUO</b>	For research use only
	Storage temperature	<b>LOT</b>	Lot
	Expiry date	<b>REF</b>	Catalogue number
	Consult instructions for use	<b>GTIN</b>	Global Trade Item Number

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